

Experiment HM-2: Electromyogram (EMG) Activity in Antagonistic Muscles

Background

The movement of parts of the body is accomplished through a system of levers composed of skeletal muscles and bones. In a lever, the muscle attached to the bone provides the effort or force that moves the bone. As the muscle contracts and relaxes, the bone, functioning as the actual lever, rotates around a joint in the skeletal system. In relation to the muscle, the bone, and the body part being moved, the joint is the fixed point that functions as the fulcrum for the lever. The body part being moved is the load in the lever (Figure HM-2-1 on page HM-2-1).

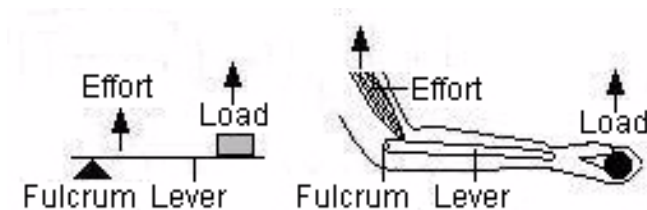


Figure HM-2-1: A Third Class lever and a counterpart in the human body. The insertion of the muscle on the bone in the forearm places the effort between the elbow (fulcrum) and the hand (load).

All levers, including the ones in the body, can be categorized into one of three classes, which are based on the position of the fulcrum in relation to the positions of the effort and the load:

- In First Class levers, the fulcrum is between the effort and the load, like on a see-saw. In the body, an example of this class is the extension of the forearm by the triceps muscle.
- In Second Class levers, the load is between the effort and the fulcrum, like a wheel barrow. In the body, an example of this class is standing on tiptoe by using the gastrocnemius muscle.
- In Third Class levers, the effort is between the fulcrum and the load, like tweezers. In the body, an example is the flexion of the forearm by the biceps muscle.

Body parts are moved in different directions by muscles that act on the same bone from different directions. A simple example is the movement of the forearm. When the biceps muscle contracts, the forearm and the hand move toward the shoulder in a motion called flexion. When the triceps muscle contracts, the forearm and hand move away from the shoulder in a motion called extension. Since these two muscles move the forearm in different directions, they are antagonistic muscles. These muscles relax and contract in a coordinated manner to place the forearm and hand in the desired position. If the effort exerted on the bone by each muscle is equal in magnitude and opposite in

direction, the forearm remains stationary. Antagonistic muscles do not need to belong to the same class of lever to work in opposition. For example, the biceps and triceps are different classes of levers; the biceps is part of a third class lever, and the triceps is part of a first class lever.

By recording the EMG activity in a muscle during the movement or positioning of a body part, it can be determined if the muscle is involved. In this experiment, students will record EMG activity from muscles on the anterior and posterior sides of the forearm to determine which ones are responsible for flexion and extension of the hand. Recording of EMG activity from these muscles will also be done while a weight is lifted by the hand. In another exercise, students will record EMG activity from the anterior and posterior sides of the lower leg to determine which muscles are active during movements or positioning that are more complex, like leaning forward, standing on toes, or rocking on heels.

Equipment Required

- PC Computer
- IWX/214 data acquisition unit
- USB cable
- IWX/214 power supply
- C-AAMI-504 ECG cable and electrode lead wires
- Disposable electrodes
- Alcohol swabs
- Small weight

IWX/214 Setup

- 1 Place the IWX/214 on the bench, close to the computer.
- 2 Check Figure T-1-1 in the Tutorial Chapter for the location of the USB port and the power socket on the IWX/214.
- 3 Check Figure T-1-2 in the Tutorial Chapter for a picture of the IWX/214 power supply.
- 4 Use the USB cable to connect the computer to the USB port on the rear panel of the IWX/214.
- 5 Plug the power supply for the IWX/214 into the electrical outlet. Insert the plug on the end of the power supply cable into the socket on the rear of the IWX/214. Use the power switch to turn on the unit. Confirm that the power light is on.

Start the Software

- 1 Click on the [LabScribe](#) shortcut on the computer's desktop to open the program. If a shortcut is not available, click on the **Windows Start menu**, move the cursor to **All Programs** and then to the listing for **iWorx**. Select **LabScribe** from the **iWorx submenu**. The [LabScribe Main window](#) will appear as the program opens.

- 1 On the **Main window**, pull down the **Settings menu** and select **Load Group**.
- 2 Locate the folder that contains the settings group, **IPLMv4.iwxgrp**. Select this group and click **Open**.
- 3 Pull down the **Settings menu** again. Select the **AntagonisticMuscles-LS2** settings file.
- 4 After a short time, **LabScribe** will appear on the computer screen as configured by the **AntagonisticMuscles-LS2** settings.
- 5 For your information, the settings used to configure the **LabScribe** software and the IWX/214 unit for this experiment are listed in Table HM-2-1 on page HM-2-2. These settings are programmed on the Preferences Dialog window which can be viewed by selecting Preferences from the Edit menu on the **LabScribe** Main window.

Table HM-2-1: Settings on the Channel Window of the Preferences Dialog Used to Configure the iWorx Recording System for Experiment HM-2.

Parameter	Units/Title	Setting	Mode/Function
Acquisition Mode		Chart	
Start		User	
Stop		User	
Display Time	Sec	10	
Speed	Samples/Sec	1000	
Channel A1	EMG Anterior	✓	3-10KHz
Channel A2	EMG Posterior	✓	3-10KHz

EMG Cable Setup

- 1 Locate the C-AAMI-504 EMG cable and electrode lead wires (Figure HM-2-2 on page HM-2-2) in the iWorx kit.



Figure HM-2-2: The C-AAMI-504 EMG cable with five lead wires attached.

- 2 Insert the black AAMI connector on the end of the EMG cable into the isolated inputs of Channels 1 and 2 of the IWX/214 (Figure HM-2-3 on page HM-2-2).
- 3 Locate the muscles of the forearm over which the recording electrodes will be placed. Muscles can be located by flexing or extending the hand and noting the areas of the forearm where the muscles are tense during these hand positions:

- One pair of recording electrodes will be placed over the flexor muscles on the anterior surface of the forearm. The first electrode in this pair will be placed about 8 centimeters below the inside of the elbow and about 4 centimeters from the medial margin of the forearm. The second electrode in this pair will be placed about 8 centimeters below the other electrode along the midline of the anterior surface of the forearm.
- A second pair of electrodes will be placed over the extensor muscles on the posterior surface of the forearm. The first electrode in this pair will be placed about 9 centimeters below the tip of the elbow along the midline of the posterior surface of the forearm. The second electrode in this pair will be placed about 6 centimeters below the first electrode and about 3 centimeters from the lateral margin of the posterior surface of the forearm.
- A fifth electrode, used as the ground, is centered between the positions of the four recording electrodes.



Figure HM-2-3: A five lead EMG cable connected to an IWX/214.

- 4 Use an alcohol swab to clean and scrub the areas where the electrodes will be placed (Figure HM-2-4 on page HM-2-2). Let the areas dry before attaching the electrodes.

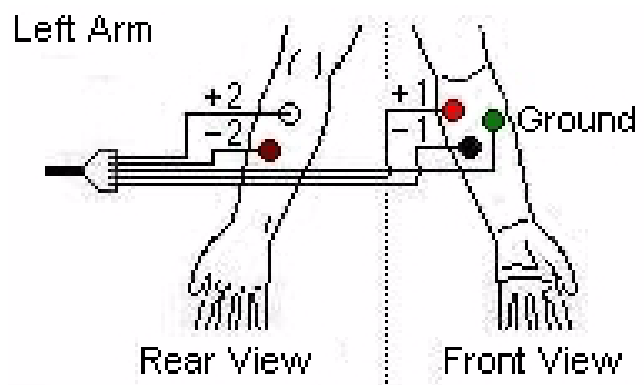


Figure HM-2-4: Position of electrodes used to record EMG from muscles in the forearm during flexion and extension.

- 5 Remove the plastic disk from a disposable electrode and apply it to one of the scrubbed areas. Attach an electrode to each of the other areas.

- Snap the recording lead wires onto the electrodes, so that:
 - the red "+1" lead is attached to the electrode on the anterior forearm that is nearest the elbow.
 - the black "-1" lead is attached to the electrode on the anterior forearm closest to the middle of the forearm.
 - the white "+2" lead is attached to the electrode on the posterior forearm that is nearest the elbow.
 - the brown "-2" lead is attached to the electrode on the posterior forearm closest to the middle of the forearm.
 - the green "C" lead (the ground) is attached to the electrode in the center of the box formed by the positions of the four recording electrodes.

Exercise 1: Antagonistic Muscles in Forearm

Aim: To study the EMG activity in muscles that work in opposition to each other to flex or extend the hand.

Procedure

- Instruct the subject that he or she will be doing the following during this exercise:
 - Before the recording begins, the subject extends his or her arm in front of their body with the palm facing upward. This position is defined as the **neutral position**.
 - Keep his or her hand open during the recording.
 - Move his or her hand upward (**flexion**) from the neutral position as far as possible and hold it in this position for four seconds (Figure HM-2-5 on page HM-2-3).

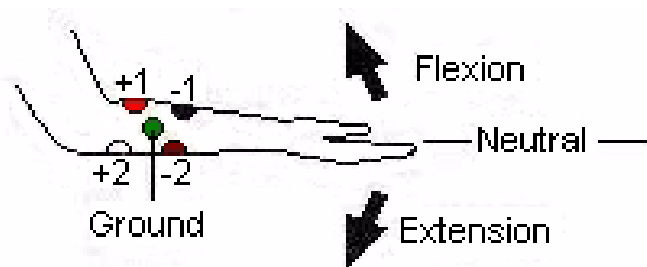


Figure HM-2-5: Movements performed while recording from muscles of the forearm responsible for flexion and extension of the hand.

- Return the hand to the neutral position for four seconds.
 - Move the hand downward (**extension**) as far as possible and hold it in this position for four seconds.
 - Return the hand to the neutral position for at least two seconds.
 - Repeat this cycle of flexion and extension two more times while recording.
- Before starting the recording, type **Neutral** in the **Mark box** to the right of the **Mark button**. Instruct the subject to place his or her hand in the **neutral position**.
 - Click the **Record button** in the upper right corner of the **LabScribe Main window**. Either click on the **Mark button** or press the **Enter key** on the keyboard to label the recording.

- While the subject's hand is in the neutral position, type **Flexion** in the **Mark box**. When the subject flexes his or her hand, click on the **Mark button** or press the **Enter key** on the keyboard to label the recording.
- While the subject's hand is in the flexed position, type **Neutral** in the **Mark box**. When the subject returns his or her hand to the neutral position, click on the **Mark button** or press the **Enter key** on the keyboard to label the recording.
- While the subject's hand is in the neutral position, type **Extension** in the **Mark box**. When the subject extends his or her hand, click on the **Mark button** or press the **Enter key** on the keyboard to label the recording.
- Repeat Steps 2 through 6 two more times.
- Repeat Steps 2 through 6 three times while the subject has formed a fist.
- When the last cycle is completed, click the **Stop button**.
- Click **AutoScale** on all channels to amplify signals. The recording should be similar to Figure HM-2-6 on page HM-2-3.
- Select **Save As** in the **File menu**, type a name for the file. Choose a destination on the computer in which to save the file, like your lab group folder. Designate the file type as ***.iwxdata**. Click on the **Save button** to save the data file.

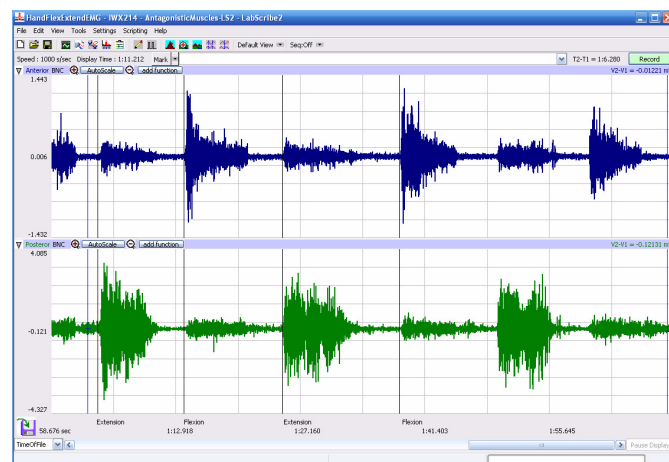


Figure HM-2-6: Recordings from anterior (top) and posterior (bottom) muscles of the forearm during flexion and extension of the hand while the subject's fingers are extended.

Data Analysis

- Scroll through the recording and find the section of data recorded while the subject was flexing and extending his or her hand.
- Use the **Display Time** icons to adjust the **Display Time** of the **Main window** so all three flexion-extension cycles appear on the **Main window**. The three flexion-extension cycles can also be selected by:
 - Placing the cursors on either side of the three adjacent cycles; and

- Clicking the **Zoom between Cursors** button on the **LabScribe** toolbar to expand the segment with the three cycles to the width of the **Main window**.

3 Click on the **Analysis window icon** in the **LabScribe toolbar** (Figure HM-2-7 on page HM-2-4) or select **Analysis** from the **Windows menu** to transfer the data displayed in the **Main window** to the **Analysis window** (Figure HM-2-8 on page HM-2-4).

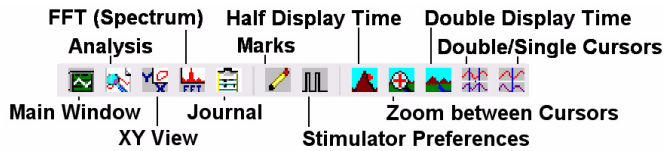


Figure HM-2-7: The **LabScribe** toolbar.

4 Look at the **Function Table** that is above the uppermost channel displayed in the **Analysis window**. The mathematical functions, **Abs. Area**, **Max-Min**, and **T2-T1** should appear in this table. The values for **Abs. Area**, **Max-Min**, and **T2-T1** on each channel are seen in the table across the top margin of each channel.

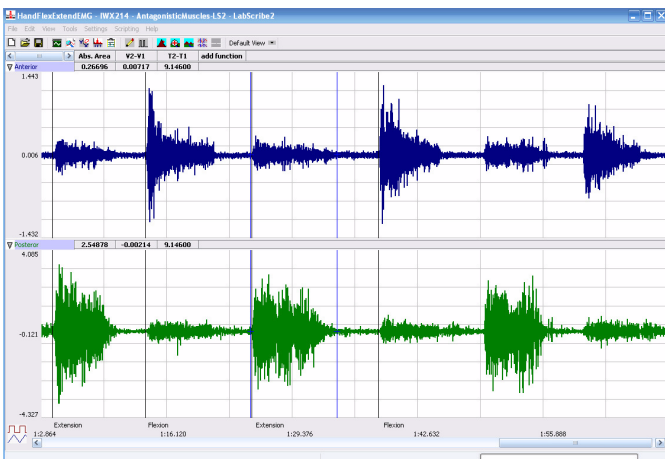


Figure HM-2-8: EMG activity from opposing muscles. Cursors are placed to measure the EMG activity taking place simultaneously in flexors and extensors.

5 Once the cursors are placed in the correct positions for measuring the three parameters of an EMG burst, the values for the parameters can be recorded in the on-line notebook of **LabScribe** by typing the names and values directly into the **Journal**, or on a data table.

6 The functions in the **channel pull-down menus** of the **Analysis window** can also be used to enter the names and values of the parameters to the **Journal**. To use these functions:

- Place the cursors at the locations used to measure the absolute areas.
- Transfer the name of the mathematical function used to determine the absolute areas to the **Journal** using the **Add Title to Journal** function in the **Anterior Channel pull-down menu**.
- Transfer the values for the absolute areas to the **Journal** using the **Add All Data to Journal** function in the **Anterior Channel pull-down menu**.

7 Use the mouse to click on and drag the cursors to the onset and offset of the EMG burst during the first flexion (Figure HM-2-8 on page HM-2-4). The values for **Abs. Area**, **Max-Min**, and **T2-T1** on the **Anterior channel** are the relative amounts of the EMG activity, the differences between maximum and minimum EMG amplitudes, and the durations of the EMG burst in the anterior muscle group during the flexion. The values for these parameters, listed in the table above the **Posterior channel**, are the same properties from the posterior muscle group during the same flexion.

8 Record the values from each muscle group during the first flexion in the **Journal**, or on a data table

9 Use the mouse to move the cursors to onset and offset of the next EMG burst, which is the first burst that occurs during extension. Measure and record the values for **Abs. Area**, **Max-Min**, and **T2-T1** from both groups of muscles during this extension.

10 Measure the **Abs.Area**, **Max-Min**, and **T2-T1** for the EMG bursts from the anterior and posterior muscles from the remaining flexions and extensions.

11 Average the values for each parameter taken from the anterior muscles during flexion. Find the average for each parameter from the anterior muscles during extension. Also, find the averages for the parameters from the posterior muscles during flexion and during extension. Enter the averages for each parameter, from each muscle group during flexion and during extension, on Table HM-2-2 on page HM-2-5.

12 Select **Save** from the **File menu**.

Questions

- Which muscles, anterior or posterior, had the most EMG activity during flexion?
- Which muscles, anterior or posterior, had the most EMG activity during extension?
- Does flexion or extension of the fingers affect the strength of EMG activity in either group of muscles?

Exercise 2: Antagonistic Muscles Doing Work

Aim: To study the EMG activity in muscles that lift weight by flexion or by extension.

Procedure

- Use the same experimental setup used in Exercise 1.
- Instruct the subject to rest his or her forearm, with the electrodes, on a flat surface with the palm up.
- Place a weight (2-3kg) in the palm of the subject's hand (Figure HM-2-9 on page HM-2-5).
- Click **Record**. Record the EMG activity from the muscles of the forearm as the subject lifts the weight by the flexion. The subject should raise and hold the weight up for four seconds and then return the weight to the table top for four seconds. Repeat this cycle two more times.
- Click **Stop** to halt the recording.

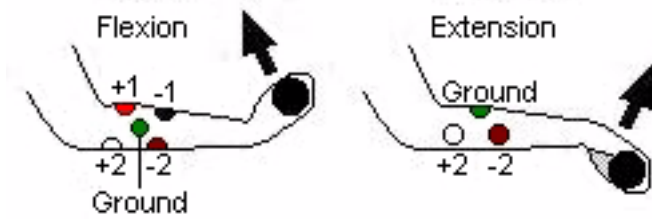


Figure HM-2-9: Hand in the palm-up position for lifting the weight by flexion (on the left), and the hand in the palm-down position for lifting the weight by extension (on the right).

- 6 Instruct the subject to rest his or her forearm on a flat surface with the palm down (Figure HM-2-9 on page HM-2-5).
- 7 Have the subject grip the same weight (2-3kg).
- 8 Click **Record**. Record the EMG activity from the muscles of the forearm as the subject lifts the weight by extension. The subject should raise and hold the weight up for four seconds and then return the weight to the table top for four seconds. Repeat this cycle two more times.
- 9 Click **Stop** to halt the recording.
- 10 Select **Save** in the **File** menu.

Data Analysis

- 1 Scroll through the recording and find the section of data recorded while the subject was lifting a weight by flexion and extension.
- 2 Use the same procedures used in Exercise 1 to measure and record the **Abs. Area**, **Max-Min**, and **T2-T1** from each muscle group while the weight was lifted by flexion and by extension.
- 3 Average the values for each parameter taken from the anterior muscles while weight was lifted by flexion. Find the average for each parameter from the anterior muscles while weight was lifted by flexion. Also, find the averages for the parameters from the posterior muscles while weight was lifted by flexion and by extension. Enter the means for each parameter, from each muscle group during flexion and during extension, on Table HM-2-2 on page HM-2-5.
- 4 Select **Save** from the **File** menu.

Questions

- 1 Does the strength of the EMG activity in the muscles of the anterior forearm differ between flexion with a weight and without a weight?
- 2 Does the strength of the EMG activity in the muscles of the posterior forearm differ between extension with a weight and without a weight?

Table HM-2-2: Relative EMG Activity in the Muscles of the Forearm during Flexion and Extension.

Forearm Action	Averages from Anterior Muscles			Averages from Posterior Muscles		
	Abs.Area	V2-V1 (V)	T2-T1 (sec)	Abs.Area	V2-V1 (V)	T2-T1 (sec)
Flexion, Hand Open						
Extension, Hand Open						
Flexion, Hand Closed						
Extension, Hand Closed						
Flexion, Lifting Weight						
Extension, Lifting Weight						

Exercise 3: Antagonistic Muscles in Lower Leg

Aim: To study muscles working in opposition to each other to maintain balance while standing.

Procedure

- 1 Locate the muscles of the lower leg over which the recording electrodes will be attached. Muscles can be located by performing dorsiflexion (lifting the toes towards the shin), or plantar flexion (pointing the toes toward the ground), and noting the areas of the lower leg where the muscles are tense during each flexion:
 - One pair of recording electrodes will be placed over the anterior tibialis muscle. This muscle is located just lateral to the tibia (shinbone) in the upper part of the calf (Figure HM-2-10 on page HM-2-6). To locate the tibialis anterior, feel for the subject's tibia. Place your fingers 2 cm to the lateral side of the

margin of the tibia and 8-12 cm below the kneecap. As the subject points his or her foot inward (supination) and upward (dorsiflexion) their foot, you should be able to see and feel the contraction of the anterior tibialis muscle beneath the skin. The first electrode in this pair will be placed about 10 centimeters below the knee on the midline of the muscle. The second electrode in this pair will be placed on the midline of the muscle, about 10 centimeters below the first electrode.

- A second pair of recording electrodes will be placed over the gastrocnemius (calf) muscle on the back of the lower leg. The first electrode in this pair will be placed about 8 centimeters below the back of the knee, along the midline of the calf muscle. The second electrode in this pair will be placed in the middle of the calf muscle along its midline, about 10 centimeters below the first electrode.
- A fifth electrode, used as the ground, is placed on the inside of the lower leg just above the ankle.

- 2 Use an alcohol swab to clean and scrub the areas where the electrodes will be placed (Figure HM-2-10 on page HM-2-6). Let the areas dry before attaching the electrodes.
- 3 Remove the plastic disk from a disposable electrode and apply it to one of the scrubbed areas.
- 4 Snap the recording lead wires onto the electrodes, so that:
 - the red “+1” lead is placed on the electrode in the upper portion of the tibialis anterior.
 - the black “-1” lead is placed on the electrode in the middle portion of the tibialis anterior.
 - the white “+2” lead is attached to the electrode near the back of the knee.
 - the brown “-2” lead is attached to the electrode in the middle of the calf muscle.
 - the green “C” lead (the ground) is attached to the electrode above the ankle.
- 5 While the subject is sitting, click **Record**. Have the subject alternate between plantar flexion and dorsiflexion of their foot. **Click AutoScale** on all channels to amplify signals. Use the techniques explained in Exercise 1 to mark the recording with comments that identify the EMG activity associated with plantar flexion and dorsiflexion.
- 6 Have the subject stand erect. Click **Record**. Have the subject rock on their feet from heels to toes and back to heels, 4 times. Mark the recording with comments that identify the EMG activity that took place.
- 7 Determine which muscle group, gastrocnemius or anterior tibialis muscles, contract when the subject rocks forward or rocks backward.

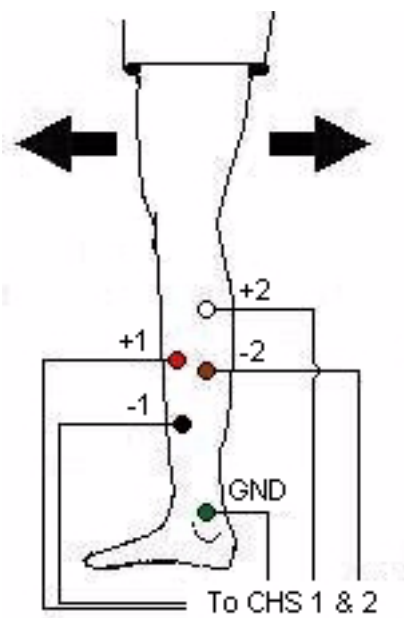


Figure HM-2-10: Position of electrodes used to record EMG from muscles in the lower leg during plantar flexion and dorsiflexion.

- 8 Have the subject rock backward and forward, 4 times, while putting more weight on the leg with the electrodes. Even though one muscle may dominate the record, the other muscle is also somewhat active. Mark the recording with comments that identify the EMG activity that took place.
- 9 Have the subject stand on one foot and remain motionless. Co-contraction of the antagonistic muscles mechanically stabilizes the joints when the subject is motionless. The stretch reflexes prevent twisting and slipping and help to maintain balance. Mark the recording with comments that identify the EMG activity that took place.
- 10 Joint stabilization is particularly important to leg and postural muscles involved in bipedal locomotion. Within the limits of artifacts induced by leg movement and cable lengths, explore the activity of the ankle flexors. Have the subject either squat or stretch upward on his/her toes. Mark the recording with comments that identify the EMG activity that took place. Click **Stop** to halt recording.
- 11 Select **Save** in the **File** menu.

Data Analysis

- 1 Scroll to the data where EMG activity was recorded from the muscles of the anterior and posterior lower leg.
- 2 Use the same procedures used in Exercise 1 to measure and record the absolute area, difference in amplitudes, and durations of the EMG bursts from the anterior and posterior muscles of the lower leg while the subject moved his or her legs and body in different positions:
 - Standing erect on both legs;
 - Rocking back and forth on both legs;
 - Rocking back and forth on the leg with the electrodes;
 - Standing erect on the leg with the electrodes;
 - Ankle flexing by squatting or standing on the toes
- 3 Enter the means for each parameter, from each muscle group during the various motions, on Table HM-2-3 on page HM-2-7

Questions

- 1 Compare the parameters for the EMG bursts from the tibialis anterior during each of the activities. When was tibialis anterior activity the greatest? The least?
- 2 Compare the parameters for the EMG bursts from the gastrocnemius during each of the activities. When was gastrocnemius activity the greatest? The least?
- 3 How does EMG activity in the gastrocnemius correlate to EMG activity in the tibialis anterior?

Table HM-2-3: Relative EMG Activity in the Muscles of the Lower Leg during Changes in Posture.

Leg Activity	Averages from Anterior Tibialis			Averages from Gastrocnemius		
	Abs.Area	V2-V1 (V)	T2-T1 (sec)	Abs.Area	V2-V1 (V)	T2-T1 (sec)
Standing Erect, Both Legs						
Rocking Forward, Both Legs						
Rocking Backward, Both Legs						
Rocking Forward, One Leg						
Rocking Backward, One Leg						
Standing Erect, One Leg						
Ankle Flex, Squat						
Ankle Flex, Standing on Toes						