


RPC-100 Respiration/Photosynthesis Chamber

Technical Note

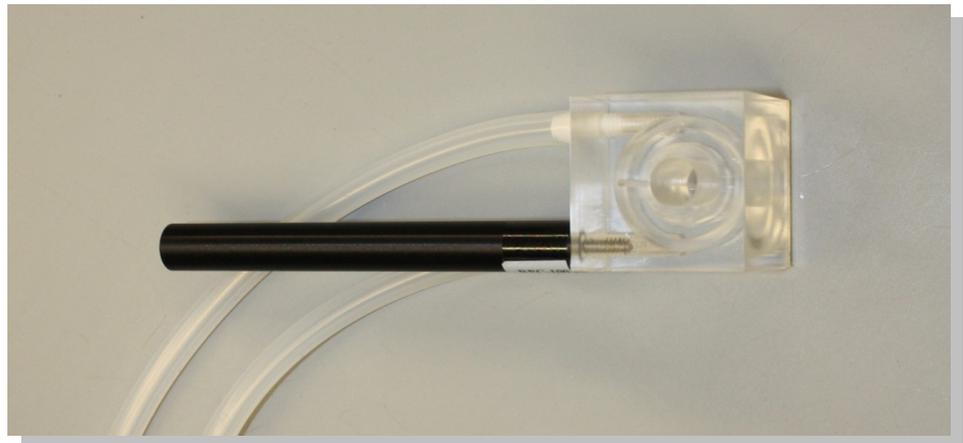

RPC-100

Overview

The RPC-100 Respiration/Photosynthesis Chamber is a reaction chamber designed for studying cellular respiration or photosynthesis. In these types of experiments, the chamber is filled with a mixture of cells or cell fractions, buffers, and reagents that react to either consume or produce oxygen. The amounts and rates of oxygen, that are consumed or produced, are measured using a dissolved oxygen electrode placed in the chamber. When mitochondria are studied, the changes in the concentration of oxygen in the chamber provide information about the movement of electrons through the chain of molecules that are responsible for the synthesis of ATP. In the typical experiments used to study mitochondria, oxygen is usually consumed. When chloroplasts are studied, the changes in oxygen concentration provide information about the light-driven movement of electrons through the chain of molecules that are responsible for creating the energy used in the synthesis of sugars. Since the experiments with chloroplasts can be performed on either their intact or their isolated photosystems, oxygen can either be produced or consumed.

How It Works

The cells or cell fractions, buffers, and reagents used in these experiments can be added or removed from the 1.6 milliliter reaction chamber of the RPC-100 through the large electrode port on the top of the chamber. The electrode port is designed to hold the iWorx ISE-730 Oxygen Electrode at the proper level in the reaction chamber during the experiment. When the electrode is removed from the chamber, the electrode port can also be used for cleaning and rinsing the reaction chamber between experiments.

The RPC-100 has a reagent port that connects to the reaction chamber from the top of the chamber. This port is smaller in diameter to restrict the flow of oxygen from the environment into the chamber. In some experiments, reagents need to be added to the reaction mixture; this can be done through the reagent port using a micropipetter with a long, thin tip.



iWorx Systems, Inc.

www.iworx.com

Labscribe is a trademark of
iWorx Systems, Inc.
©2015 iWorx Systems, Inc.

RPC-100 Respiration/Photosynthesis Chamber

During experiments, the reaction mixture needs to be mixed to promote the exchange of oxygen across the membrane of the electrode. This is accomplished by placing a small magnetic stir bar in the bottom of the reaction chamber. This stir bar is controlled by a stir motor placed under the chamber.

Since some experiments need to be conducted at temperatures other than room temperature, the temperature of the reaction mixture can be maintained by circulating heated or cooled water through the jacket that surrounds the reaction chamber.

How to Use the RPC-100

Equipment Setup

- 1) Attach the RPC-100 chamber to a ring stand or similar support device.
- 2) Place a stir motor under the chamber.
- 3) If the temperature of the reaction is going to be maintained at a temperature other than room temperature, connect the tubing between the water jacket and the optional circulating pump.
- 4) Rinse the reaction chamber several times with deionized water before calibrating the oxygen electrode.
- 5) Fill the reaction chamber with room temperature deionized water in preparation for the calibration of the oxygen electrode.
- 6) Place a small, clean Teflon stir bar in the bottom of the reaction chamber and align the stir motor to place the stir bar in the center of the reaction chamber.
- 7) Rinse the membrane and membrane holder of the ISE-730 Oxygen Electrode several times with deionized water before putting it in the electrode port.
- 8) Carefully place the ISE-730 in the electrode port of the chamber.

Experiments

Labscribe experiments using the RPC-100 include:

- **Experiment CM-5: Carbon Dioxide Fixation in Intact Cells** (found in the **Cellular Metabolism** category of the Labscribe **Settings** menu as **CO2Fixation**)
- **Experiment CM-3: Mitochondrial Respiration** (found in the **Cellular Metabolism** category of the Labscribe **Settings** menu as **MitoRespiration**)
- **Experiment CM-4: Photosynthesis in Isolated Thylakoids** (found in the **Cellular Metabolism** category of the Labscribe **Settings** menu as **Photosynthesis**)



iWorx Systems, Inc.

www.iworx.com

RPC-100 Respiration/Photosynthesis Chamber

Technical Data and Specifications

SPECIFICATIONS	
Construction	Clear acrylic plastic
Volume of Reaction Chamber	1.6 milliliters
Designed for use with	ISE-730 Oxygen Electrode and DO2-100 Current-to-Voltage Adapter
Also Included	2mm x 7mm Teflon coated stir bars (2) Gel-loading micropipette tips



iWorx Systems, Inc.

www.iworx.com

iWorx Systems, Inc. 62 Littleworth Road, Dover, New Hampshire 03820
(T) 800-234-1757 / 603-742-2492 (F) 603-742-2455